

All data taken at Pacific Northwest National Laboratory (PNNL)
Operators: Steven W. Sharpe, Timothy J. Johnson and Robert L. Sams : sw.sharpe@pnl.gov
Version 1.0, February, 03

Composite spectrum for DMDS_50T

Effective burden of composite spectrum: 1 part-per-million-meter (ppm-meter) at 296 K

Equivalent concentration x path-length of composite spectrum: 3.8778×10^{-6} grams/liter-meter

Contaminated with 0.11% of 2,3-Butadione (out gassing from previous sample) and corrected for by spectral subtraction of 2,3-Butadione followed by rescaling the composite spectrum.

Sample Conditions-

- Chemical name and CAS number: Dimethyldisulfide, 2,3-dithiabutane, sulfa-hitech, methyl disulfide, (methyldithio)methane, DMDS, $\text{CH}_3\text{SS}(\text{CH}_3)$: [624-92-0]
- Physical properties: fw=94.1894 g/mole, fp=-85° C, bp=109.7° C
- Supplier and stated purity: Aldrich, 99+%
- Sample class: I (PNNL scale).
- Temperature of sample: 49.94 ± 0.02 C
- Diluent: Sample back filled with ultra high purity nitrogen to 760 ± 5 Torr
- Individual samples at 1.06950, 4.0893, 16.24, 2.2995, 8.4848 and 20.28 Torr. Path length = 19.96 cm. Final data is a composite spectrum.
- Preparation: Multiple freeze-thaw cycles at -90 C to remove air.

Instrument Parameters-

- Bruker-66V FTIR, temperature controlled environment, evacuated optics bench
- Modified to include second aperture, between interferometer output and sample cell. This substantially reduces both “ghosting” and warm aperture effects.
- Spectral range: 6,500 to 600 cm^{-1} (1.534 to 16.667 microns)
- Instrumental resolution based on maximum interferometer displacement is 0.112 cm^{-1}
- Spectral interval after 2X zero-filling interferogram and FFT: 0.06 cm^{-1}
- Interferogram zero-fill: 2X
- Apodization: Boxcar
- Phase correction: Mertz
- Beam splitter: Potassium bromide (KBr)
- IR source: Carbide glowbar (22 V)
- Scanner velocity: 60KHz (HeNe crossing frequency)
- Number of interferograms averaged per single channel spectra: 256
- Detector: Mid-band HgCdTe, photoconductive, 77K operation
- Folding limits: 15798 to 0 cm^{-1}

Post Processing and Related Parameters-

- Non-linearity detector correction (Bruker proprietary) applied to interferogram ($\tau = 0.85$, $\nu = 530$)
- Composite spectrum created from 6 individual absorbance (base-10) spectra via classical least squares fit: Intercept=0, slope is fitted, individual absorbance values weighted by T^2 (transmission squared), all absorbance values > 1.6 are given zero weight
- Calculated and estimated errors: Type A = 0.69%, Type B = 3%
- Frequency correction (already applied): $V(\text{corrected}) = V(\text{instrument}) * 0.99999896 + 8.812 \times 10^{-4}$
- Axis units: X=wavenumbers (cm^{-1}), Y=Absorbance (base-10)
- Baseline correction via 7th order polynomial subtraction